# **Operating instruction**



# Top Loading Balances "L/LW" series

Filename : MAN\_L\_EN\_S3\_00\_R1\_0.doc January 2011

#### CONTENTS

1	INSTALLATION INSTRUCTION	
2	STORAGE	4
3	WEIGHING PAN ASSEMBLY	5
4	KEYPAD AND DISPLAY	6
5	OVERVIEW	7
	5.1 FRONT PART OF THE BALANCE, L / LW MODELS	
	5.2 REAR AND BOTTOM PART OF THE BALANCE, L/LW MODELS	
6		
	<ul><li>6.1 STAND BY</li><li>6.2 SIMPLE WEIGHING</li></ul>	
7		
/	7.1 External calibration	
	<ul> <li>7.1 EXTERNAL CALIBRATION</li> <li>7.2 INTERNAL CALIBRATION</li> </ul>	
	7.2.1 Autocalibration (AUT-CAL)	
	7.2.2 Internal calibration (I-CAL)	
	7.2.3 External calibration (E-CAL)	
	7.2.4 Technical calibration (TEC-CAL)	
8	TARE FUNCTION	
	8.1.1 Manual tare function (Not available for V-range models)	
	8.1.2 Optional alphanumeric external keyboard (code T201)	
9		
1(	0 PC COMMUNICATION SETTING	
11	1 PRINTER SERIAL COMMUNICATION SELECTION	
12	2 TRANSMISSION SPEED SELECTION	
13	3 AUTOZERO FUNCTION	
14	4 FILTERS SELECTION	
15	5 STABILITY FUNCTION	
16	6 CONTRAST REGULATION	
17	7 LCD BACKLIGHT SETUP	
18	8 AUTO POWER-OFF FUNCTION	
19	9 PIECECOUNTING FUNCTION	
	19.1 VISUALIZATION OF TOTAL AND UNIT WEIGHT OF PIECES	
	19.2 MANUAL INSERTION OF THE UNIT AVERAGE WEIGHT	
	19.3 AUTOMATIC UPDATE OF UNIT WEIGHT	
20	0 V-RANGE FUNCTION	
21	1 DENSITY DETERMINATION OF A SOLID OR A LIQUID	

	1.1 1.2	DENSITY DETERMINATION OF A SOLID	
22	MA	XIMUM LOAD DETERMINATION FUNCTION	9
23	PER	RCENTAGE WEIGHING FUNCTION 4	0
	3.1 3.2	Mode with weight reference	-
24	ANI	IMAL WEIGHING FUNCTION 4	3
25	TO	TALIZING 4	4
26	TH	RESHOLD FUNCTION 4	6
2	6.1 6.2 6.3	BOTH THRESHOLDS SET	7
27	RS2	32 INTERFACE FEATURES 4	8
28	WO	DRKING WITH BATTERY*	3
29	ERF	ROR CODES	4
30	MA	INTENANCE AND CARE	5
31	QUI	ICK GUIDE TO BALANCE PARAMETERS SETUP5	6
32	QUI	ICK GUIDE TO THE USE OF THE BALANCE PROGRAMS5	7
<b>33</b> DE	<b>BAI</b> FINIT	L <b>ANCE TECHNICAL CHARACTERISTICS</b> ERRORE. IL SEGNALIBRO NON 1 10.	È
34	ACC	CESSORIES	8
35	WA	RRANTY	9
36	EQU	UIPMENT DISPOSAL	9
37	COI	NFORMITY DECLARATIONERRORE. IL SEGNALIBRO NON È DEFINITO	).


# **1** Installation instruction




#### WARNING:


Please follow carefully these steps for installing and use the new balance before starting your work routine. A way of use of the instrument different from this user manual will not guarantee the instrument's safety anymore.


- **Remove** the balance, calibration weight( if included) and accessories from the cartoon and check for any visible damage of the instrument.
- **Do not install** the balance in a place with air flows, heavy thermal changes and vibrations.
- Do not use the balance in blast risk environment.
- The humidity rate of the balance environment must be between 45% and 75%

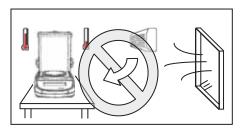

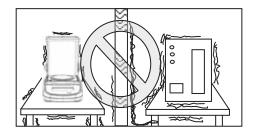

- **Place** the weighing pan and the support pan on the balance (see par. 3).
- Level the balance using the level bubble and levelling feet located underneath the case

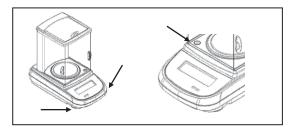

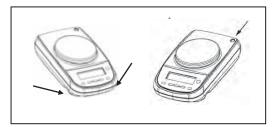


- Connect power supply to connector 2 located on the rear panel of the unit (see par. 5).
- Insert power unit into AC outlet, which shall be easily accessible; after few seconds the balance will automatically switch on.
- Wait 30 minutes from switch on and then calibrate the balance using the appropriate mass, following the instructions (par. 7)


- Calibrate again the balance every time it's moved to another place
- **Check** balance calibration periodically.
- We recommend not dropping heavy objects on balance pan, in order to avoid damages.
- Assistance service must be effected by specialized staff and the spare parts used must be original. Therefore, it is necessary to apply to the reseller who sold the equipment.

#### 2 Storage


- Storage temperature: +5 °C...+40°C
- Storage humidity 45% 75%.
- Keep package in the case the balance should be sent back to the factory for service. Disconnect the cables and the accessories for avoiding damages during transportation.
- **Do not place** the balance in extreme temperature and humidity condition and avoid the balance to take strong hit.

# 3 Weighing pan assembly

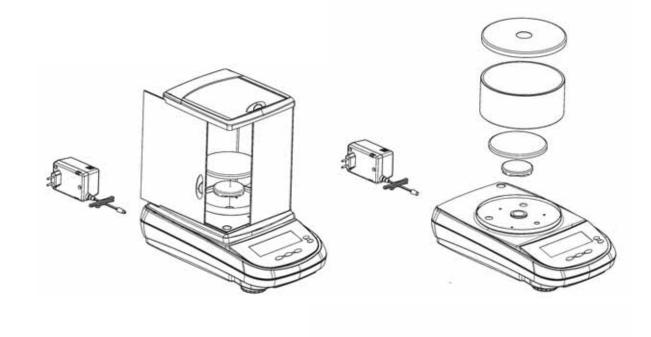

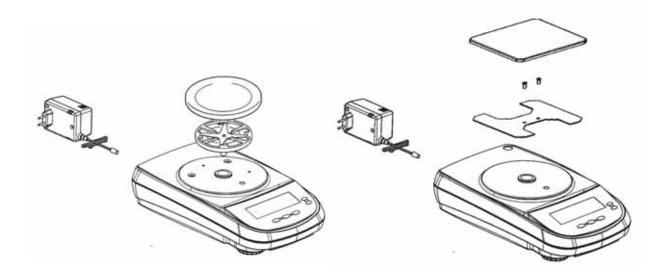


# 4 Keypad and display

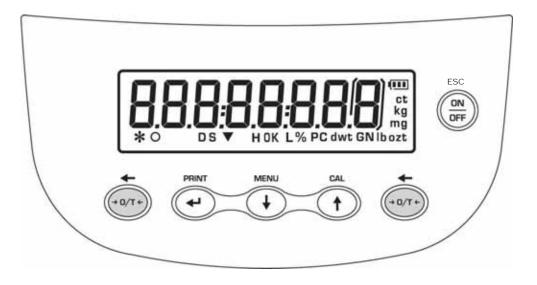

ESC ON OFF	Standby (OFF/ON) or escape ESC.	*	Stability indicator
		0	Zero indicator
(→ 0/T ←	TARE or zero button.	%	Percentage weighing
PRINT		PC	Piece counting
(+)	Selection CONFIRM or SEND data to printer.		Battery charge indicator
	·		Insert data mode
MENU	Balance setup MENU button.	н	Upper threshold
		L	Lower threshold
CAL	Balance CALIBRATION button.	DS	Density measure mode
		ct, ozt, l	Measure unit <b>b, GN, dwt, Kg, mg</b>

# 5 Overview

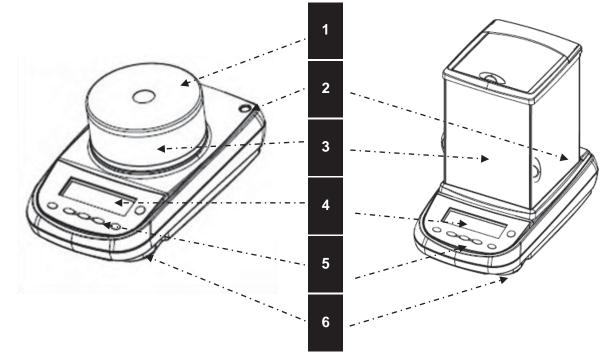


### 5.1 Front part of the balance, L / LW models

Fig.1 Balance with circular and three-panes windshield

Nar	me and function
1	Cap of windshield
2	Levelling bubble
3	Windshield
4	LCD display
5	Function buttons
6	Levelling feet

### 5.2 Rear and bottom part of the balance, L / LW models

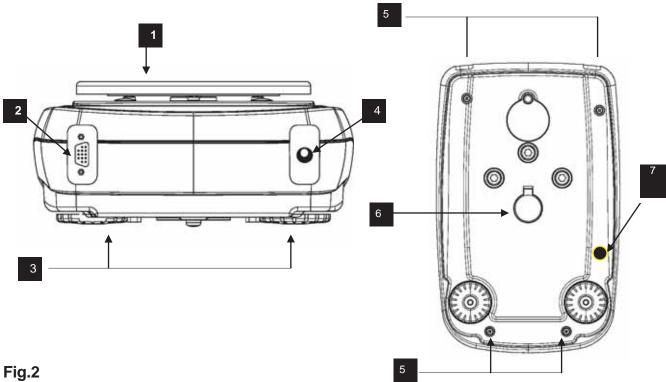


Pannello posteriore e parte inferiore bilancia

Nar	me and function
1	Weighing pan
2	9 poles (pin) female connector for RS232 interface for printer/PC/external keyboard
3	Levelling feet
4	connector for power supply
5	Screws for balance closing. In models with 4 feet it is necessary first to remove the
5	two rear feet (one fixed and one moving) to access to the closing screws
6	Hook for under-balance weighing. Hook not available in internal calibration models.
Ŭ	Note: please remove the cap to access the hook
	Protection screw for autocalibration motor (only for models with internal calibration).
7	ATTENTION: remove this screw by hand as soon as the balance is removed from
	its package

# 6 Weighing

After having connected the balance to AC outlet, it will perform an internal circuits test, after that the balance will set itself in stand-by mode.

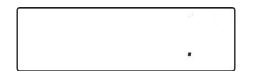

#### 6.1 Stand By


From "STAND BY" mode:

- Press **ON/OFF** button to bring balance to work conditions.
- Press again ON/OFF button to return to "STAND BY" condition.



#### 6.2 Simple weighing

Load the sample to weigh on the pan and read the value on display as soon as the stability symbol  $\mathbf{x}$  (*star*) appears

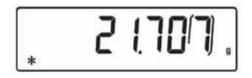

# 7 Calibration

Electronic balances take mass measurements making use of gravity (g). Difference of latitude in geographic areas and altitude will vary gravity acceleration value (g). Therefore, for accurate measurements, the balance must be adjusted to the local environment. This adjustment is accomplished by calibration function.


#### 7.1 External calibration

Calibration is accomplished by pressing **CAL** button.

1. Press **CAL** button when pan is empty, dashes are displayed on the display.




2. When calibration weight value starts to flash, load the weight on the pan.

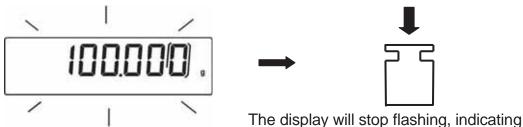

calibration weight value.

Once the calibration is effected will be shown the value of the calibrated weight and the current unit of measure.


 Unload calibration weight from the pan. The balance is ready for weighing operations.

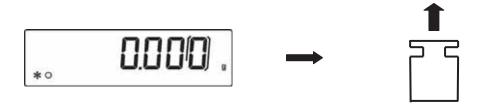

NOTE: if there is interference during calibration process, an error message will be displayed.

Moreover, it is possible to calibrate the balance with a calibration weight higher than the one set by default:

a) Press and keep **CAL** button pressed with empty pan until the acoustic alarm stops, then release the button. On display it will be visualized the string "-**CAL**-", followed by flashing string "LOAD".

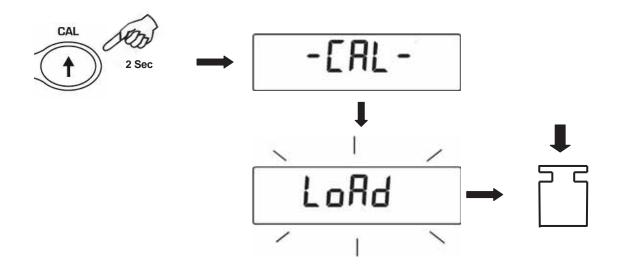

b) Load on the pan a weight equal higher or lower than default calibration weight; the balance will recognize it as valid weight if equal or higher than calibration weight as long as it is a whole number in comparison with the most meaningful digit of calibration weight.

e.g.: if calibration weight is 200g, it will be possible to calibrate the balance with values from 100g 200g, 300g, 400g up to the highest limit of balance weighing range. The message "LOAD" on display will stop flashing. Once calibration has been effected, the value of calibrated weight will be displayed.

c) Unload calibration weight. The balance is ready for weighing operations.



NOTE: if there is interference during calibration process, an error message will be displayed.


#### 7.2 Internal calibration

\* only for models with internal calibration "-i"

In these balance models there are 4 calibration modes:

From display zero condition, press and keep pressed the **MENU** button until the acoustic alarm gets mute, then release the button. The message "**unitS**" will be visualized on display, press then **MENU** button until you visualize "**Calib**" on display. Press **PRINT** to confirm.

- 1. Select the calibration mode you wish by pressing **MENU** button in sequence:
  - AUT-CAL: auto calibration
  - I-CAL: internal calibration
  - E-CAL: external calibration
  - TEC-CAL: technical calibration

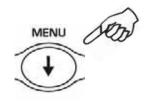

- Press PRINT button to confirm "AUT-CAL", "I-CAL", "E-CAL". To confirm "TEC-CAL" keep pressed the PRINT button until the acoustic alarm gets mute.
- 3. After selection, the balance returns to calibration menu. Press and keep pressed **MENU** button until the acoustic alarm gets mute, then release the button. Balance is again ready for weighing operations.

#### 7.2.1 Autocalibration (AUT-CAL)

The balance autocalibrates through the internal reference calibration mass, after the microprocessor has checked that no other weighing operations are being effected. In this mode it is furtherly possible to effect calibration through reference internal mass by pressing **CAL** button at any time, being sure first that there's no weight load on the pan.

 Press CAL button with empty pan. The message "CAL" will be displayed and balance calibration will be effected automatically.



2. At the end of calibration, balance returns to normal weighing conditions.



If calibration is not finished due to vibrations or air flows, then the message "CAL

**bUt**" is displayed. Press again **CAL** button. If the problem repeats then select external calibration and contact the product supplier.

#### 7.2.2 Internal calibration (I-CAL)

The balance calibrates itself with internal reference mass <u>ONLY</u> on user command by pressing **CAL** button.

Before effecting internal calibration, be sure that there's no weight loaded on the balance pan.

#### 7.2.3 External calibration (E-CAL)

The balance is calibrated through an external reference mass (**NOT** provided with models with internal calibration).

(Follow procedures described at paragraph 7.1)


#### 7.2.4 Technical calibration (TEC-CAL)

This function allows storing the value of internal reference mass whenever checking or assistance actions require it.

1. After having selected the **TEC-CAL** calibration mode, press **CAL** button at empty pan. It will be displayed "CAL".



2. When the value of calibration weight start flashing on display, load the weight on to the balance pan.

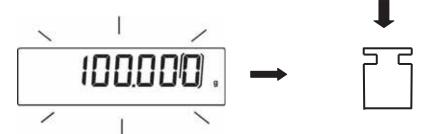

- 3. Wait the acoustic alarm and that the displayed calibrated weight value stops flashing, then unload the weight from balance pan.
- 4. When string "**0.000**" is displayed continuously, then press and keep pressed the **PRINT** button. This starts the internal weight value automatic acquisition and store. During the acquisition cycle, the display will show "**TEC-MEM**".




- 5. After having stored the value of internal calibration weight, balance returns to normal weighing conditions.
- 6. Return to calibration menu as described at paragraph 6.2 and set the desired calibration mode: internal, automatic or external..



<u>ATTENTION</u> : this procedure must be effected only using E2-class reference masses.

### 8 Tare function

1. Load the container on the pan. The display will show the weight.

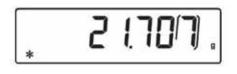



2. Press O/T button. "O-t" string will be displayed



3. After reaching stability, the value "**0.000**" will be displayed. If the stability is not reached (due to air flows or vibrations or other disturbs) the dashes will remain displayed.



4. Load the objects to weigh in the container. Read net weight value on display.






#### 8.1.1 Manual tare function (Not available for V-range models)

This function allows to insert manually the tare values.

- 1. Press and keep pressed the **O/T** key whit no objects on the weighing pan until the beep alarm stops, then release the key.
- 2. Will be shown on the display the following string:

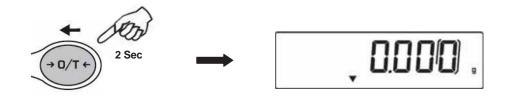


3. Insert now the desired tare value using the keys **CAL** and **MENU** for increase or decrease the value, and press the **O/T** key for skip to the following digit. During the value insert mode, if keep pressed the **O/T** key is possible to delete the inserted value.



4. Press **PRINT** key to confirm the value inserted.


#### 8.1.2 Optional alphanumeric external keyboard (code T201)

If you have the optional external alphanumeric keyboard, then it is possible to perform tare also by pressing the **TARE/DELETE** button of this keyboard, in the same way as previously described.

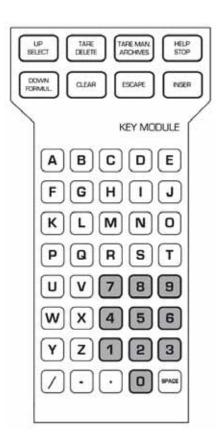

It's also possible to manually insert a known tare value by the keypad.

1. Press **TARE MAN** button on the alphanumeric keyboard. An arrow will appear on the balance display along with the previous value of manual tare, if one has been inserted before.

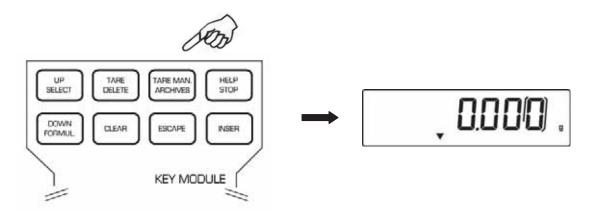

2. Press **CLEAR** to set at zero the previous value (if any) then insert the new value using the alphanumeric buttons placed on the bottom part of alphanumeric keyboard.


- 3. Press **INSER** to confirm.
- 4. Press **ESCAPE** button to escape from tare condition.

To set to zero the value of the manual tare inserted by means of alphanumeric keyboard, effect a normal operation of tare by pressing **O/T** button of the balance or with **TARE/DELETE** button placed on the optional alphanumeric keyboard.


# 9 Weight units

It is possible to select the weight unit that balance will use to display weight.

1. From display zero condition press and keep pressed the **MENU** button until the acoustic alarm gets mute, then release the button. The message "**unitS**" will be displayed then press **PRINT** to confirm.

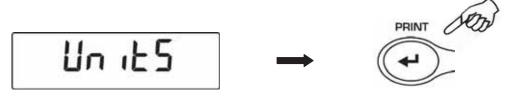


2. It will be displayed "**GRAM**" unit. Pressing now the **MENU** or **CAL** button, it will be possible to scroll forward or backward the weight units menu.

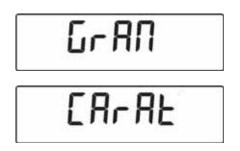


3. Press **PRINT** button to confirm or **MENU** button to shift to the other weight unit.

	SYMBOL	UNIT	CONVERSION FACTOR 1g =
NOW	GRAM	GRAM	1.
CAL COO	CARAT	CARAT	5.
$( \mathbf{A} )$	OUNCE	ONCE	0.035273962
	POUND	POUND	0.0022046226
	PENN.	PENNYWEIGHTS	0.643014931
	ONCETR.	ONCE TROY	0.032150747
	GRANO	GRAIN	15.43235835
- I	TAEL HON	HONG KONG TAEL	0.02671725
MENU COD	* TAEL SGP	SYNGAPORE TAEL	0.02646063
	* TAEL ROC	R.O.C. TAEL	0.02666666
$\checkmark$	* MOMME	MOMME	0.2667
	M 10	x 10	10.
ŧ	M 100	x 100	100.

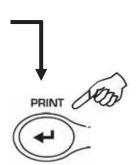

- 4. After having selected the desired weight unit, press the **MENU** button to go to next parameter or **CAL** button to return to previous.
- 5. To escape from parameters setup menu, press the **MENU** button until the acoustic alarm gets mute, then release the button. The balance returns to normal weighing conditions.


\* The unit of measure with the star (\*) are not available in the balances with the software for paper grammage. In these models is possible to set from the menu of the units of measure the multiply factor of the weight "M 10" or "M 100". For visualizing the m<sup>2</sup> weight, select the M 100 multiply factor (as described at par.9) and load on the pan a 10x10 cm piece of paper.

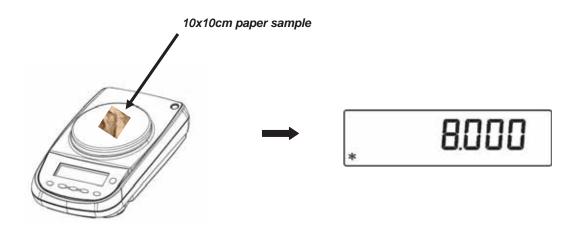


# **10 PC communication setting**

Connect the balance to PC with proper cable (par. 25)

 From zero display condition press and keep pressed the MENU button until the acoustic alarm gets mute, then release the button. It will be displayed the string "unitS", then press the MENU button until it is displayed "PC-PRTR" message then confirm by pressing PRINT button.

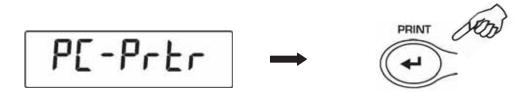

 Press the MENU button until the "PC cont" is displayed to choose continuous print or "PC CMd" to choose print to PC at user command, then press the PRINT button to confirm the choice.

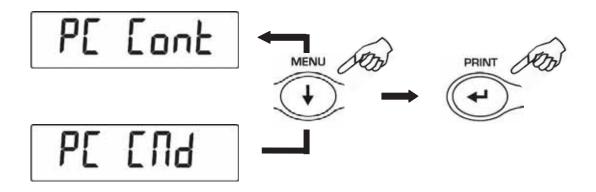

- 3. After having selected the desired transmission mode, press the **MENU** button to go to next parameter or the **CAL** button to go to previous.
- 4. To escape from parameters setup menu, press the **MENU** button until the acoustic alarm gets mute, then release the button. The balance returns to normal weighing conditions.
- 5. The balance returns to normal weighing conditions sending the data continuously or at user command.

*0	

NOTE: select speed transmission (par. 12)


### **11** Printer serial communication selection

Connect the balance to the printer using the proper cable (par. 25)


 From zero condition on display, press and keep pressed the MENU button until the acoustic alarm gets mute, then release the button. The message "unitS" will be displayed, then press MENU until you read the message "PC-PRTR" on display and confirm pressing PRINT button



- To select the data print mode, press MENU until the message "PRINT" is displayed or "TLP50" if you want to connect the TLP50 printer for printing on labels or on paper with date and time.
- 3. Press **PRINT** to confirm.

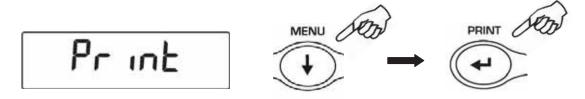

- 4. After having selected the data print mode, press **MENU** button to go to next parameter or **CAL** button to go to previous one.
- 5. To escape from parameters setup menu, press **MENU** button until the acoustic alarm gets mute, then release the button. The balance returns to normal weighing conditions, ready to transmit data each time the **PRINT** button is pressed.

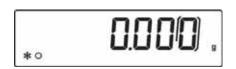


NOTE: select speed transmission (par. 12)

# 12 Transmission speed selection

1. From zero condition on display, press and keep pressed the **MENU** button until the acoustic alarm gets mute, then release the button. The message "**unitS**" is displayed, then press **MENU** button until the message "**BAUD RT**" is displayed and confirm by pressing the **PRINT** button.



2. Select serial data transmission speed (1200-2400-4800-9600 baud). Pressing **MENU** or **CAL** buttons it will be possible to scroll forward or backward trough the different transmission speeds, then confirm your choice by pressing **PRINT** button.

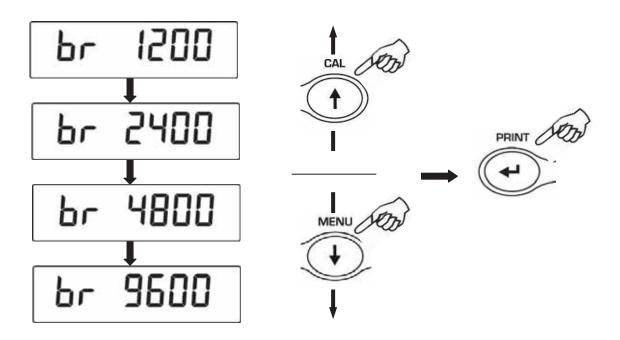

- 3. After having selected the transmission speed you wish, press the **MENU** button to go to next parameter or **CAL** button to go to previous one.
- 4. To escape from parameters setup menu, press the **MENU** button until the acoustic alarm gets mute, then release the button.
- 5. The balance will return to standard weighing conditions.




# **13 Autozero function**

Autozero is an automatic correction of a possible zero drift.

- Au0 OFF = autozero disabled
- Au0 1 = soft autozero
- Au0 2 = medium autozero
- Au0 3 = heavy autozero
- Au0 3E = heavy autozero over all range
- From zero condition on display, press and keep pressed the MENU button until the acoustic alarm gets mute, then release the button. The message "unitS" is displayed, then press MENU button until the message "AUTO 0" is displayed, then press PRINT to confirm.



 Pressing MENU or CAL button it will be possible to scroll forward or backward through the different autozero levels, select the one you wish and confirm it by pressing the PRINT button.

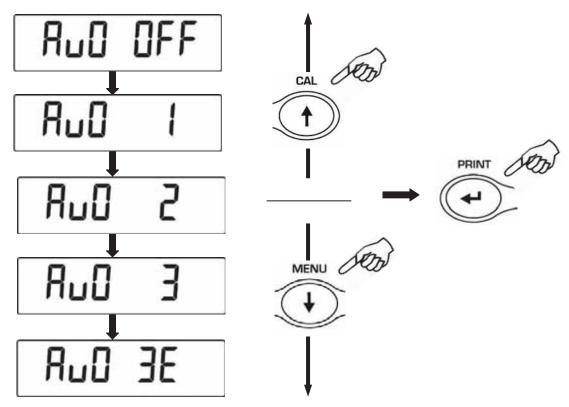

- 3. After having selected the autozero level you wish, press the **MENU** button to go to next parameter or **CAL** button to go to previous one.
- 4. To escape from parameters setup menu, press the **MENU** button until the acoustic alarm gets mute, then release the button.
- 5. The balance will return to standard weighing conditions.


# 14 Filters selection

It is possible to adapt the balance to the different environment conditions thanks to the selection of three filters:

- FILTER 1: proportion of ingredients condition
- FILTER 2: stable conditions
- FILTER 3: unstable conditions
- 1. From zero condition on display, press and keep pressed the **MENU** button until the acoustic alarm gets mute, then release the button. The message "**unitS**" is displayed, then press **MENU** button until the message "**FILTER**" is displayed then confirm it by pressing the **PRINT** button.



2. Pressing **MENU** or **CAL** button it will be possible to scroll forward or backward the different filtering levels, select the one you wish and then confirm it by pressing the **PRINT** button

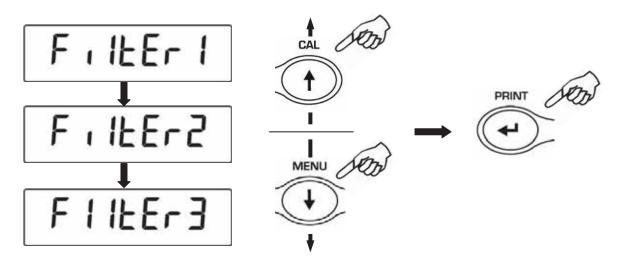

- 3. After having selected the filtering level you wish, press the **MENU** button to go to next parameter or **CAL** to go to previous one.
- 4. To escape from parameters setup menu, press the **MENU** button until the acoustic alarm is over, then release the button.
- 5. The balance will return to standard weighing conditions.




NOTE: It is suggested to use FILTER 1 when proportion of ingredients must be performed

# **15 Stability function**

The stability symbol is displayed when the weight is stable inside a defined range

- **STAB 1** = for stable environments
- STAB 2 = for not so stable environments
- **STAB 3** = for unstable environments
- 1. From zero condition on display, press and keep pressed the **MENU** button until the acoustic alarm gets mute, then release the button. The message "**unitS**" is displayed, then press **MENU** button until the message "**StAbiL**" is displayed, then confirm this by pressing the **PRINT** button.
- 2. Pressing **MENU** or **CAL** button it will be possible to scroll forward or backward the different stability levels, select the one you wish and then confirm it by pressing the **PRINT** button.

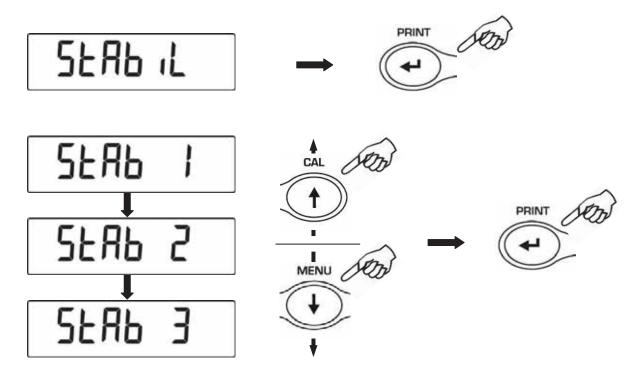

- 3. After having selected the stability level you wish, press the **MENU** button to go to next parameter or the **CAL** button to go to previous one.
- 4. To escape from parameters setup menu, press the **MENU** button until the acoustic alarm gets mute, then release the button.
- 5. The balance returns to normal weighing conditions.



# **16 Display Contrast regulation**


It is possible to regulate the contrast level of display to have a comfortable view of the indication at different angles of usage.

There are 15 different levels of regulation:

 From zero condition on display, press and keep pressed the MENU button until the acoustic signal gets mute, then release the button. The message "unitS" is visualized, now press the MENU button repeatedly until the message "contr" is displayed and confirm this by pressing the PRINT button



2. Now, pressing the buttons **MENU** or **CAL** it is possible to increment or decrement the level of the display contrast; select the one desired and confirm by pressing the **PRINT** button.

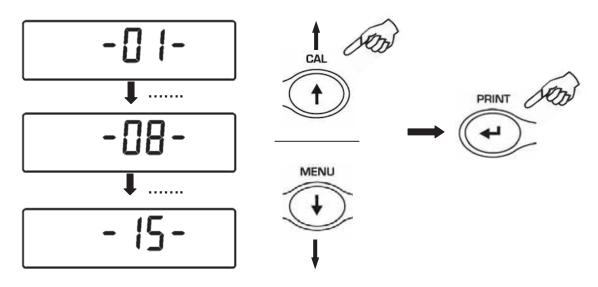

- 3. After you have confirmed the contrast level desired, press the **MENU** button to go to the next parameter or the **CAL** button to go to previous one.
- 4. To escape from parameters setup menu, press the **MENU** button until the acoustic alarm gets mute, then release the button.
- 5. The balance returns to normal weighing mode



# 17 Display Backlight setup


The balance display is equipped with backlight to make the indication more visible also during low light conditions.

There are 3 working modes:

- **ON** = light always switched ON
- **OFF** = backlight always switched OFF
- AUTO = backlight automatically switched on during weighing operations
- From zero condition on display, press and keep pressed the MENU button until the acoustic alarm gets mute, then release the button. The message "unitS" is displayed, then press MENU button until the message "bLt", then press the PRINT button to confirm this.



2. Pressing **MENU** or **CAL** button it will be possible to scroll forward or backward the different working modes, select the one you wish and then confirm it by pressing the **PRINT** button

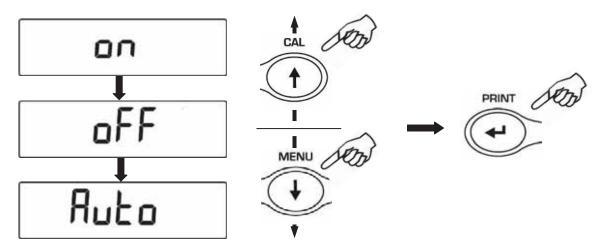

- 3. After having selected the backlight working mode you wish, press the **MENU** button to go to next parameter or the **CAL** button to go to previous one.
- 4. To escape from parameters setup menu, press the **MENU** button until the acoustic alarm gets mute, then release the button.
- 5. The balance returns to normal weighing conditions.



# **18 Auto Power-off function**

This function allows to activate the automatic power-off of the balance after a defined balance idle-time. There are 4 auto power-off modes:


- disab = Auto power-off disabled
- **2 Min** = Auto power-off after 2 minutes of idle time

#### 5 Min = Auto power-off after 5 minutes of idle time

- 15 Min = Auto power-off after 15 minutes of idle time
- From zero condition on display, press and keep pressed the MENU button until the acoustic alarm gets mute, then release the button. The message "unitS" is displayed, then press MENU button until the message "tiME oFF" then confirm by pressing the PRINT button.



2. Pressing **MENU** or **CAL** button it will be possible to scroll forward or backward the different auto power-off modes, select the one you wish and then confirm it by pressing the **PRINT** button

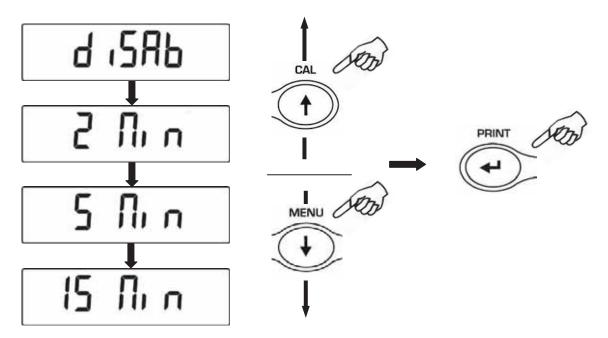

- 3. After having selected the auto power-off mode you wish, press the **MENU** button to go to next parameter or the **CAL** button to go to previous one.
- 4. To escape from parameters setup menu, press the **MENU** button until the acoustic alarm gets mute, then release the button.
- 5. The balance returns to normal weighing conditions.

### **19** Piececounting function

1. From zero condition on display, press **MENU** button until the message "**Count** " is displayed then press the **PRINT** button to confirm



Select the number of pieces to put on the pan as sample by pressing MENU repeatedly to increase and CAL to decrease the number. Choice of the number of pieces (10, 25, 50, 100) depends from the weight of single pieces. If available, load the empty container first then press PRINT to confirm.



Load on the pan the number of pieces displayed on the balance display

LoAd



Press again the **PRINT** button and wait the weight to stabilize.


If the number of samples is enough (for example 10 as in figure), then this number will be displayed and it will be possible to go on by loading the pieces to count on the balance pan.



If the weight of the pieces to count is too low in comparison with balance resolution, then an error message will be displayed. In this case it is necessary to use a balance with higher resolution. If the weight of samples is acceptable but not enough, then the message "**Add**" will be displayed.

Add a quantity of pieces so that to almost double the number read on display, then press the **PRINT** button

If the number of pieces is still not enough then the message "Add SMP" is displayed. Add again a quantity of pieces so that to almost double the number you read on display

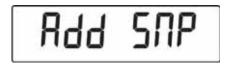




As soon as it is reached a sufficient number of pieces, this number is displayed and it will be possible to go on with count by loading the pieces count on the balance pan.



3. To escape from piececounting mode press the **ON/OFF** button and the balance returns to normal weighing conditions.

#### 19.1 Visualization of total and unit weight of pieces

1. Press **MENU** button to pass from pieces number visualization to total weight visualization.

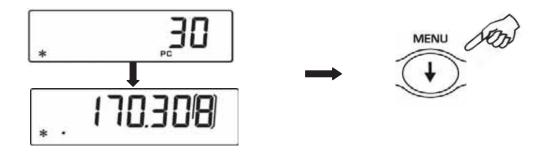

- 2. In order to display the weight of single piece from the total number of pieces press and keep pressed the **MENU** button until the acoustic alarm gets mute.
- 3. Press again the **MENU** button to display the number of pieces.

#### 19.2 Manual insertion of the unit average weight

It is possible to insert the unit average weight of the sample, if known. In this way can be avoided the sampling of the units.

 From zero condition on display, press the MENU button until the message"Count", then press the PRINT button to confirm. Select MANUAL by pressing repeatedly the MENU button.



2. Press the **PRINT** button to confirm.



3. Insert the unit weight of a sample in grams using **CAL** and **MENU** buttons for increase and decrease the value and use the **O/T** button to skip to the others digits. For insert the For inserting the decimal point keep pressed the **CAL** button. Keeping pressed the O/T button is possible to delete the inserted value.



- Press the **PRINT** button to confirm.
   If the inserted weight is 100 times lower than balance resolution, it will be displayed an error message.
   To escape without inserting the weight press **ON/OFF.**
- 5. If the weight is enough then "**0**" is displayed ; it is now possible to go on to count loading the pieces on the pan.
- 6. Press **MENU** button to display the total number of pieces, then press again the button to display again the number of pieces.
- 7. To escape from piececounting function, press **ON/OFF** button.

It is possible to use the optional alphanumeric keyboard to insert the weight of the unit sample as following:

 From zero condition on display, press the MENU button until the message"Count", then press the PRINT button to confirm. Select MANUAL by pressing repeatedly the MENU button.



- 2. Press the **PRINT** button to confirm.
- Insert the unit weight of a sample in grams using the numerical buttons from 0 to 9 and the decimal point.
   In case of mistake, press the CLEAR button to start insertion again.
- 4. Press the **INSER** button to confirm.
- If the inserted weight is 100 times lower than balance resolution, it will be displayed an error message.
   To escape without inserting the weight press ESCAPE (on alphanumeric keyboard) or ON/OFF.
- 6. If the weight is enough then "**0**" is displayed ; it is now possible to go on to count loading the pieces on the pan.
- 7. Press **MENU** button to display the total number of pieces, then press again the button to display again the number of pieces.
- 8. To escape from piececounting function, press **ON/OFF** button.

#### 19.3 Automatic update of unit weight

After having effected the sampling, it is possible to update the average piece weight as follows:

- 1. Instead of loading all the pieces to count, load a number of pieces equal to the double of the ones loaded on the pan and wait for the acoustic alarm.
- It is now possible to repeat this procedure up to a maximum of 255 pieces or to proceed to the normal counting of the pieces. This procedure allows a more accurate evaluation of the average unit weight and a better precision in the pieces count.


NOTE: automatic update is not active if the sampling has been effected by insertion of the average unit weight

# 20 V-Range function

\* This (optional) function is available ONLY for double range models \*

 V-RANGE function is to be considered when it is necessary to effect proportioning of small quantities with accuracy. To activate V-RANGE function, after having put on the pan the container if any, press keep pressed the TARE button until the acoustic alarm gets mute.

In this way a tare operation is automatically performed, at the end of which it is possible to operate with the highest resolution range.

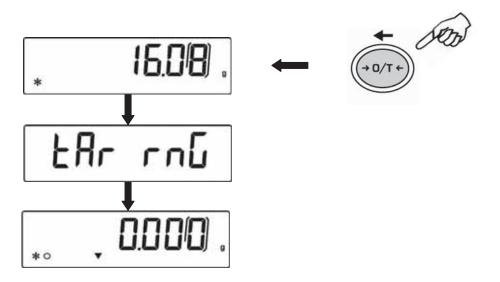

- 2. The range change is pointed out by the switch-on of a small arrow on the display.
- The range of highest sensivity is kept active as far as the net weight exceeds in positive the maximum value set for the lower sensitivity range. However it is possible to disable this function pressing and keeping pressed the TARE button until the acoustic alarm gets mute. In both the cases total weight is displayed.
- 4. In V-Range function the manual tare insertion by (optional) alphanumeric keyboard is not available.

# 21 Density determination of a solid or a liquid

\* The balance must be equipped with the proper optional software (not available for all models) \*


The software for density determination allows to calculate the density of a solid or a liquid through the under balance hook or through the optional hydrostatic kit (cod. T027)

#### 21.1 Density determination of a solid

1. From zero condition on display, press the **MENU** button until the message "**dEnS**" is displayed, then press the **PRINT** button to confirm.



2. Select the function "d SoLid" through the MENU button and then confirm with **PRINT** button.

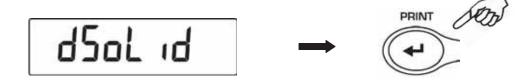

3. The density of the liquid to use will be displayed, the default value is equal to 1.0000 (distilled water at 20°C).



 It is possible to set a different value using the CAL e MENU buttons for increase and decrease the value and using the O/T button for skip to the others digits. Keeping presses the O/T button is possible to delete the inserted value.



- 5. It is also possible to set a different value if the balance is equipped with the optional alphanumeric keyboard.
- 6. Once the wished value is set, press **PRINT** button.
- 7. Now it will be asked to weigh the solid in air, confirm this with **PRINT** button.




- 8. If necessary, effect a tare operation and load the solid. Press the **PRINT** button to acquire the value. During value acquisition the message **WEI AIR** will be flashing.
- Now it is asked to weigh the solid inside the liquid. Effect the tare of the small basket inside liquid, immerse the solid and press **PRINT.** During the acquisition of the value the message **WEI LIq** will be flashing.



10. The result of density of the solid is then displayed. Pressing the **PRINT** button it is possible to print the value of density, if the balance is equipped with a printer.



11. If any error occurs, this message will be displayed:

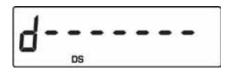

12. Press now the **ON/OFF** button to escape from density function or the **MENU** button to perform another measurement.

### 21.2 Density determination of a liquid

1. From zero condition on display, press the **MENU** button until the message "**dEnS**" is displayed, then press the **PRINT** button to confirm.



2. Using the MENU button, select the function "d Liquid" and confirm this with PRINT



3. It will be displayed the value of the density of the glass float of known volume to use, the default value is equal to 3.0000.




 It is possible to set a different value using the CAL e MENU buttons for increase and decrease the value and using the O/T button for skip to the others digits. Keeping pressed the O/T button is possible to delete the inserted value.



- 5. It is possible to set a different value if you have the optional alphanumeric keyboard
- 6. Once the wished value is set, press the **PRINT** button.

7. It is now asked to weigh the glass float in air, confirm this with **PRINT** 

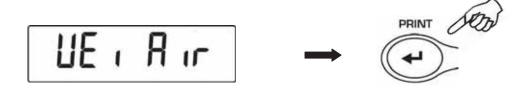

- 8. Now perform a tare if necessary and put the glass float to the weighing hook. Press then the **PRINT** button to acquire the value. During the acquisition of the value the message **WEI AIR** will be flashing.
- It is now asked to weigh the glass float immersed into the liquid. Perform a tare if necessary and immerse the glass float into the liquid. Press the **PRINT** button. During the acquisition of the value the message **WEI LIq** will be flashing.



10. The result of the density of the liquid is now displayed. Pressing the **PRINT** button it is possible to print the value of density, if the balance is equipped with a printer.

11. In case of error the following message is displayed

12. Press the **ON/OFF** button to escape from density function, or the **MENU** button to perform another measurement.

# 22 Maximum load determination function

\*The balance must be equipped with the specific optional software (not available for all models)\*

The function **M LOAD** allows to measure the maximum load for a solid.

1. To activate the function M LOAD from zero display condition, press repeatedly the **MENU** button until the message "**M Load**" is displayed, then press the **PRINT** button to confirm. When this function is activated, a tare is automatically performed.



2. Activation of this function is indicated by the visualization of the letter **M** on the last digit on the display.



3. It is now possible to determine the weight of the maximum load.



4. Then press the **TARE** button to effect another measurement.



- 5. When the function **M LOAD** is active, the calibration is disabled.
- 6. Press the button **ON/OFF** to escape from **M LOAD** function.

# 23 Percentage weighing function

This function allows to read the weighing value as a percentage of a reference weight. The reference weight is took as 100% percentage value (factory setting). There are two modes for acquiring the reference weight: one is automatic (with the reference weight) and one manual (by inserting manually the value of the reference weight).

### 23.1 Mode with weight reference


1. From zero condition of the display press the **MENU** button until will be displayed the "**Perc**" message, then press the **PRINT** button to confirm.



2. Press the MENU button to select the "**Perc A**" function and press **PRINT** button to confirm.




3. Tare function will be performed and will be shown on the display the string "Load". Then load on the weighing pan the reference and press the PRINT button, the string "Load" will start to flash and once the weight is acquired will be visualized the value with the % indication.

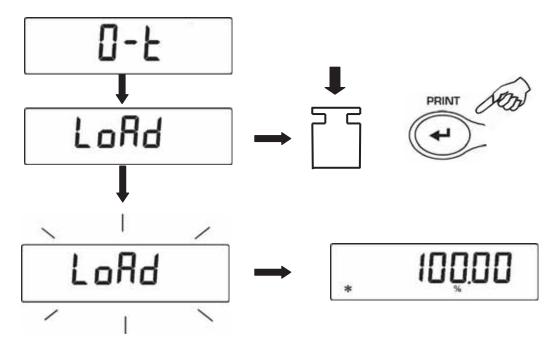

4. Remove now the reference weight and load the sample. Is possible to read the weight in percentage now.

5. Press the **MENU** button to visualize the weight in gram and vice-versa.

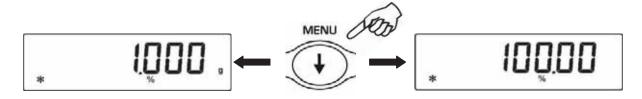

6. Press the **ON/OFF** button to escape from the percentage weight function.


NOTE: If the reference weight is lower than 10 digits after the acquiring sequence of the weight, the string ERROR 07 will be displayed.

### 23.2 Mode with manual insertion of the reference weight.

1. From zero condition of the display press the **MENU** button until will be displayed the "**Perc**" message, then press the **PRINT** button to confirm.



2. Press the MENU button to select the "**Perc A**" function and press **PRINT** button to confirm.

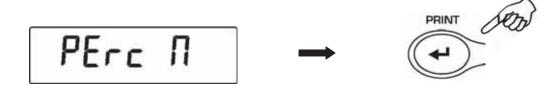


 It is possible to set a different value using the CAL e MENU buttons for increase and decrease the value and using the O/T button for skip to the others digits. Keeping pressed the O/T button is possible to delete the inserted value. The inserted value will be stored in the memory until the balance is switched off.



- 4. Press **PRINT** button when the desired value of the reference weight is inserted
- 5. Load now the sample and read the percentage value.



6. Press the **MENU** button to visualize the weight in gram and vice-versa.

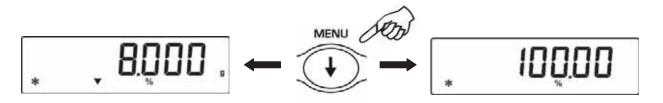


- 7. Press the **ON/OFF** button to escape from the percentage weight function.
- 8. It is also possible to set the value if the balance is equipped with the optional alphanumeric keyboard. Press PRINT button to confirm the inserted value.

NOTE: If the reference weight inserted is lower than 10 digit displayed, will be shown the ERROR 07.


# 24 Animal weighing function

This function allows to acquire the averaged weight of moving objects or animals for a time period that can be set.


1. From zero condition of the display press the **MENU** button until will be displayed the "anim UEi" message, then press the **PRINT** button to confirm.

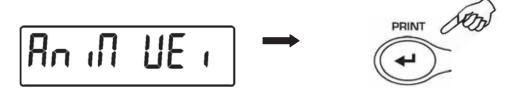

2. Set the desired period time from 5 to 90 sec. using the button **MENU** to decrement and **CAL** to increment time.Confirm by pressing the **PRINT** button.

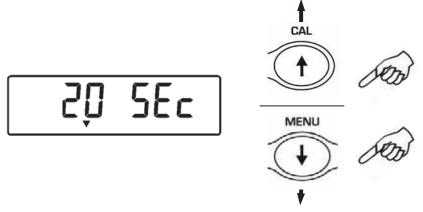

3. It is now displayed the value of current weight, effect tare if necessary, then load on the pan the sample to weight and press the **PRINT** button

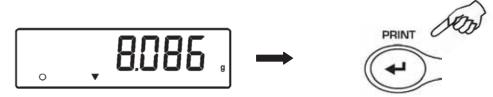

- 4. It is then displayed the count-down of the sampling time value that has been previously set.
- 5. When the count-down is finished, it is calculated and then displayed the value of the averaged weight measured.



6. Press once the **ON/OFF** button to start another measurement, or twice to escape this function and return to weighing mode


# 25 Totalizing

This function allows to effect sum of consecutive weighings.


1. From zero condition of the display press repeatedly the **MENU** button until will be displayed the "**add**" message, then press the **PRINT** button to confirm.




2. The Tare is automatically effected.

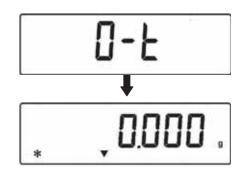

3. Load the first weight to sum.

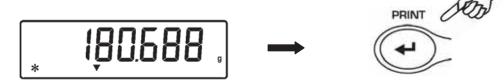


- 4. Confirm with the **PRINT** button.
- 5. The Tare is again effected.

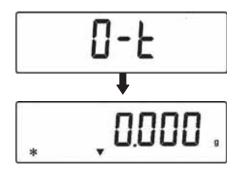


6. Load the second weight to sum.



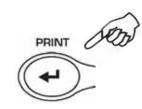

PRINT

- 7. Confirm with the **PRINT** button. Follow all the above steps for all weights to sum up to a maximum number of 99.
- 8. Pressing the **MENU** button it is possible to show in sequence the informations respectively of total weight, number of weighings effected and current weight.

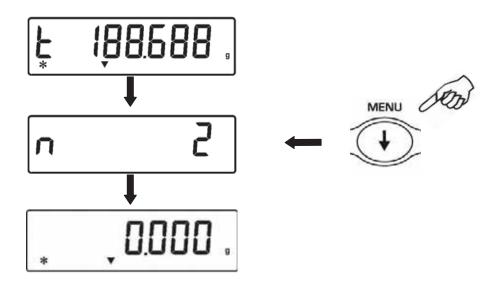

9. Press the **ON/OFF** button to escape this function and return to weighing mode

# 26 Threshold function

Threshold function allow to determinate if the weight loaded on the pan is above or below the two threshold fixed by the user.


1. From zero condition of the display press the **MENU** button until will be displayed the string "**H o L**", then press the **PRINT** button to confirm.



 Insert the lower threshold value using the CAL e MENU buttons for increase and decrease the value and using the O/T button for skip to the others digits. Keeping pressed the O/T button is possible to delete the inserted value. The inserted value will be stored in the memory until the balance is switched off.



- 3. Insert the higher threshold value as same as the lower value steps.
- 4. The string "**bEEP OFF**" will be displayed and select by the **MENU** button if the acoustic alarm should be set or not if the weight is within the two threshold values set before. Press **PRINT** button to confirm.

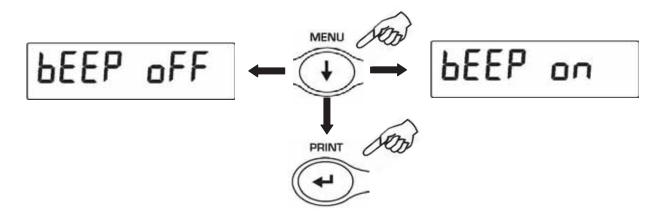

 If the value of the thresholds is correctly inserted, the balance will be back to the weighing mode with the indication of the threshold state (H = higher threshold, L = lower threshold, OK= weight within the two set thresholds).

### NOTE: If the values are not inserted correctly will be displayed the string ERROR 07.

There are 3 modes of the threshold function:

### 26.1 Both thresholds set

This mode allow to set the lower and the higher threshold values and to identify the acceptable range for the weight value shown by the switching on of the symbol "OK" and by the acoustic alarm is activated. When the value of the weight is under the lower threshold value set will be displayed the symbol "L" while if the weight is over the higher threshold value set will be displayed the symbol "H".

### 26.2 Lower threshold set only

Setting only the lower threshold and setting to zero the higher threshold, will be displayed the symbol "OK" and eventually the acoustic alarm if activated whenever the loaded weight is higher than the lower threshold value set. When the value of the weight is under the lower threshold set the symbol "L" will be displayed.

### 26.3 Higher threshold set only

Setting only the higher threshold and setting to zero the lower threshold, will be displayed the symbol "OK" and eventually the acoustic alarm if activated whenever the loaded weight is lower than the higher threshold value set. When the value of the weight is above the higher threshold set the symbol "H" will be displayed.

# 27 RS232 interface features

### 1. General features

The balance transmits the value visualized on the display following serial RS232C standard, allowing to print the value of weight to a PC monitor or to a serial printer. In the case of connection to a PC, it will be possible to select the transmission in continuous mode or transmission at user command through pressing of the **PRINT** button ( as described at par.10). The balance is also capable of receiving commands, always through the standard RS232C, that allow performing all the functions available through the keyboard of PC itself. The speed of transmission and reception can be selected, as described previously (see par.12), to 1200, 2400, 4800, e 9600 baud. The character format is of 8 bit preceded by one bit of start and followed by a bit of stop. Parity is not considered.

### 2. Selection of interface for PC

Selecting the transmission to PC (personal computer IBM compatible), it is achieved a continuous transmission output, at the same rate of weight update on display of the balance. It is possible to perform all the functions of the balances directly from the computer keyboard, transmitting to the balance the ASCII codes as in the table below. The connector to use for connection to PC is the number 2 with reference to fig.2 of paragraph 5 of this manual.


CODE	1 <sup>st</sup> FUNCTION (SINGLE PRESS)
"T" = H54	TARE
"C" = H43	CALIBRATION
"E" = H45	ENTER
"M" = H4D	MENU
"O" = H4F	ON/OFF

CODICE	2 <sup>nd</sup> FUNCTION (PROLONGED PRESS)
"t" = H74	TARE
"c" = H63	CALIBRATION
"e" = H65	ENTER
"m" = H6D	MENU
"o" = H6F	ON/OFF

Selecting the transmission to PC at user command, it is achieved a transmission output only when the **PRINT** button is pressed, also in this case it is possible to perform all the functions of the balance directly from the keyboard of the computer, sending to the balance the ASCII codes in the table above.

### 3. Connection of the balance to PC

To receive/transmit data, link the connector (n.2 of fig.2 at par.5) of the balance to the serial port of your Personal Computer as shown below:

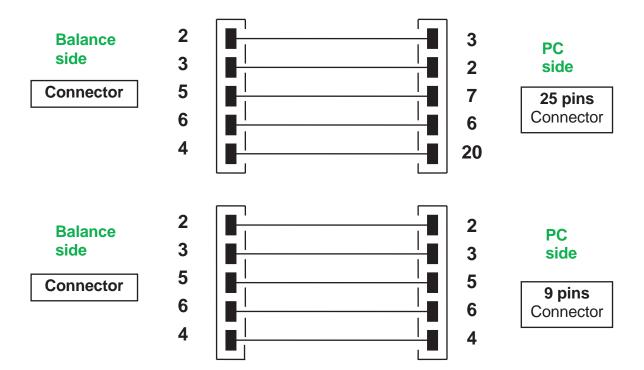

### 4. Transmission format

String transmitted is composed by the following 14 characters:

- First character: weight sign (blank or -)
- second/ninth character: weight or other data
- tenth/twelfth character: weight unit symbol
- thirteenth character: stability indicator
- fourteenth character: carriage return
- fifteenth character: line feed

Eventual non-significative zero is spaces.

In the following tables the various transmission formats are shown:

Weighing mode (valid both for continuous transmission and transmission at user command)

1°	2°	3°	4°	5°	6°	7°	8°	9°	10°	11°	12°	13°	14°	15°
Sign				We	ight				W	'eight u	nit	Stability	CR	LF

Density mode (only in transmission at user command mode)

1°	2°	3°	4°	5°	6°	7°	8°	9°	10°	11°	12°	13°	14°	15°	16°	17°
d	=		D	Dens	sity v	valu	е		Space		V	Veight uni	t		CR	LF

### Piececounting mode (only in transmission at user command mode)

### Number of pieces

1°	2°	3°	4°	5°	6°	7°	8°	9°	10°	11°	12°	13°	14°	15°	16°
	Pcs		:		spa	ces				Nur	nber	of pie	ces		

### Total weight of pieces

1°	2°	3°	4°	5°	6°	7°	8°	9°	10°	11°	12°	13°	14°	15°	16°	17°	18°	19°	20°
		We	ight			:	space			Va	alue o	f wei	ght			space	g	space	S

### Average unit weight of pieces:

	<u> </u>																	
1°	2°	3°	4°	5°	6°	7°	8°	9°	10°	11°	12°	13°	14°	15°	16°	17°	18°	
	PMU		:		spa	ces				Va	alue o	f weig	ght			spaces	g	

PMU stands for average unit weight

Percentage weight mode (only in transmission at user command mode)

### Percentage

1°	2°	3°	4°	5°	6°	7°	8°	9°	10°	11°	12°	13°	14°	15°	16°	17°	18°
	Pe	erc				space	9				Perce	entage	Ð			sapce	%

Weight

1°.	2°	3°	4°	5°	6°	7°	8°	9°	10°	11°	12°	13°	14°	15°	16°	17°	18°
		We	ight			spa	ace			V	Veigh	t valu	е			sapce	g

Animal weighing mode (only in transmission at user command mode

Time	е																
1°	2°	3°	4°	5°	6°	7°	8°	9°	10°	11°	12°	13°	14°	15°	16°	17°	18°
	Tir	ne		space	=		space	<u>.</u>	Time	value		Sec			spa	ace	

Averaged Weight

_																	
1°.	2°	3°	4°	5°	6°	7°	8°	9°	10°	11°	12°	13°	14°	15°	16°	17°	18°
	Ave			=		spa	ace		Å	Vera	ged v	veight	value	Э	spa	ace	g


### 5. Selection of interface to printer

Selecting the PRINTER mode, the serial output of the balance is set to work with serial printers.

In this case the printing is effected only when the **PRINT** button is pressed and with stable weight. If the stability is not reached within ten seconds, the message **ERR05** is displayed preceded by a short acoustic signal and the value of the weight is not sent to the printer.

### 6. Connection of the balance with the serial printer

Connect a serial printer to the connector (n.2 of fig.2 at par.5) to the balance as shown in the following scheme:

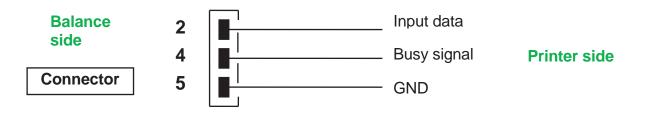

If the optional TLP50 printer is used, it will be possible to print both in *continuous-mode* and in *labels-mode* with the following formats:

Weighing mode and maximum loa	d mode
-------------------------------	--------

12-02-2008	12:00
Weight:	22.000 g

### **Piececounting mode**

12-02-20	008 12:00	
Pcs Weight: PMU:	100 300.000 g 3.000 g	

Density determination mo	de l	Percentag	e weigh	t mode
12-02-2008 12:00		12-02-20	09	12:00
d= 2.80066 g/cm3d		Perc. Weight:	100. 300.00	0% 0 g

### Animal weighing mode

Time = 6 Sec
Ave. = 59.446 g

12-02-20	09	12:00
Perc. Weight:	100 300.00	


### Totalizing (sum of weights) mode


12-02-2	2009	12:00
1. 2.  99.	16.589 17.226	g g
33.		
S=	33.815	g

### 7. Connection of the balance to the optional alphanumeric keyboard

To connect the optional external alphanumeric keyboard to the balance, it must be used the same connector that is used to link the balance to PC. In this case the connection to PC or to printer must be effected through the connector placed on the optional alphanumeric keyboard.

8. Map of RS232 interface Connector (n.2 fig.2 par.5)

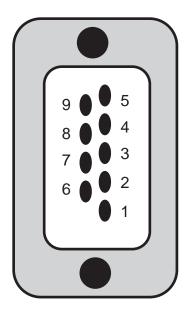

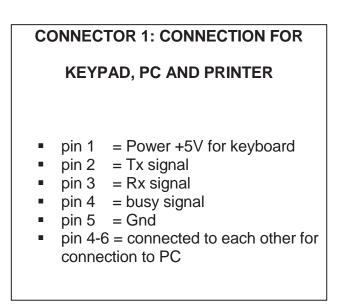




# 28 Working with battery\*

\*This operation mode is not available for the models with internal calibration

The balance can work, at wish, using only the battery power. The battery on board is located inside the balance and grants about 30 hours of autonomy

- With the external power supply disconnected, it is possible to switch on or switch off the balance trough the **ON/OFF** button and work without being connected to power AC outlet.
- 2. When the balance is working using the battery, the symbol of battery with charge level is displayed on up-right corner of the balance's display. The three lines all switched on means a maximum level of battery charge, instead when all three lines are switched off and only the battery symbol is displayed means that it is necessary to connect the balance to AC power outlet to re-charge the battery.

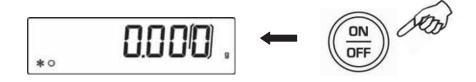

- 3. To re-charge the battery, connect the provided external power supply to the balance. (See par.5)
- 4. When the external power supply is connected, the balance automatically switch itself on and, after performing the initial test on display, it enters the stand-by mode showed by a single point in the display. In standby mode also the battery recharge symbol will remain switched on.

1	

5. The charge level will step by step raise until it reaches the maximum level. However let the battery recharge at least for 8-10 hours before disconnecting the external power supply from the balance.

NOTE: Battery duration depends also from working mode of the display backlight. To get a longer battery duration, switch off the backlight or set it in AUTO mode.

ERROR MESSAGE ON DISPLAY	MEANING	POSSIBLE SOLUTIONS
ERR01	Weight not stable after operation of tare	Protect the balance from air flows or from vibrations of the working table
ERR02	impossible to start the calibration due to instability of the balance	Protect the balance from air flows or from vibrations of the working table.
ERR03	calibration weight not correct or balance unstable	Calibrate with correct weight or protect the balance from environment disturbs
ERR04	weight of samples for the piececounting function not adequate or unstable	Select a bigger number of samples or protect the balance from vibrations.
ERR05	impossible to print / weight unstable	Protect the balance from environment disturbs
ERR06	Weight cannot get stable in density mode	Protect the balance from environment disturbs
ERR07	error in inserting the data / weight unstable	
ERR08	Anomaly on autocalibration motor	Contact service staff
"UNLOAD"	weight loaded on the pan or pan not positioned properly	Remove the weight from the pan or position properly the pan and underpan.
"CAL But":	the balance requires to be re- calibrated	Unload weights, if any, on the pan, and press the CAL button
ry	Over-range condition	Unload the weights loaded on the pan
L	Under-range condition	Place properly pan and underpan

## 30 Maintenance and care

Regular maintenance of yours balance guarantee accurate measurements.

### Cleaning

Before cleaning the balance unplug the power supply of the balance from the voltage supply of your room. Do not use aggressive cleaning product (as solvents or similar), use a humid towel with soft detergent, Avoid liquids to go inside the instrument during the cleaning. Wipe the balance with a soft towel. Parts of samples or powder can be removed using a brush or vacuum cleaner.

### Safety checks


Safety of the instrument is no more guaranteed when: -balance power supply is clearly damaged

-balance power supply is not working anymore

-balance power supply is stored for long time in hard environment conditions. In these instances refer to the assistance centre where specialized technician will make reparations to bring back the instrument in the safety conditions eventually.


31 Quick guide to balance parameters setup

- To enter the balance parameters setup menu, press and keep pressed the MENU button until the acoustic alarm gets mute.
- Use then the MENU button to go to next parameter, use the CAL button to go to previous and the PRINT button to confirm the choice.
  - To escape from menu, press and keep pressed the MENU button until the acoustic alarm gets mute.

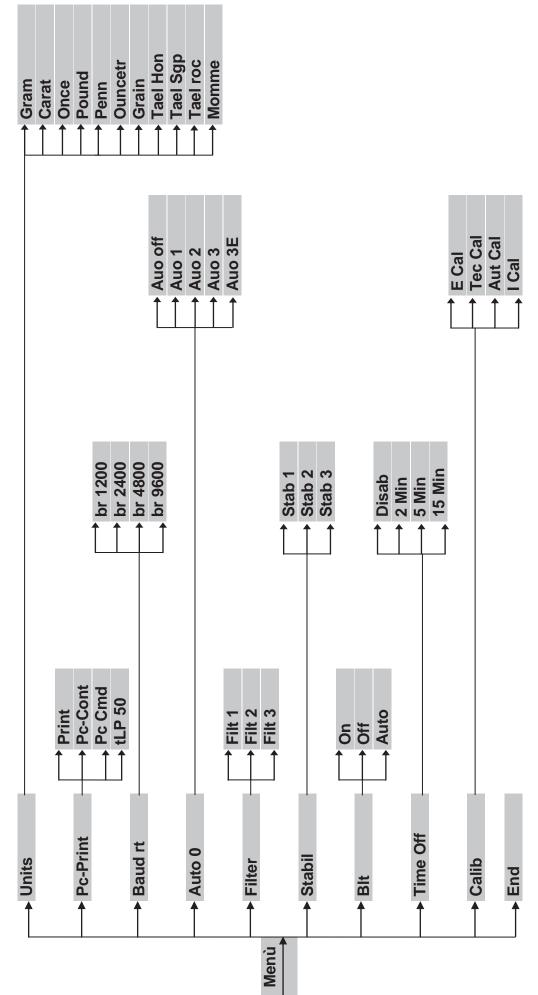



- To enter the balance programs menu, press the MENU button.
- Use then the **MENU** button to go to next parameter, use the **CAL** button to go to previous and the **PRINT** button to confirm the choice. To escape from menu, press and keep pressed the **MENU** button until the acoustic alarm gets mute.

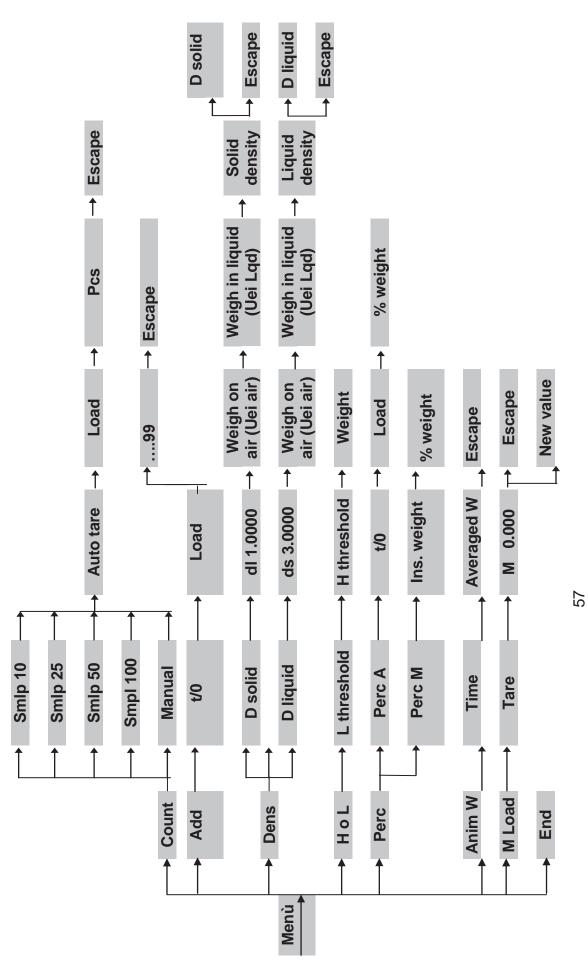

# 33 Optional accessories

ACCES	SORIES
Code	DESCRIPTION
T201	EXTERNAL KEYOBARD mod. KEY-MODULE
T002	SOLID DENSITY KIT
T027	SOLID LIQUID DENSITY KIT
T214	RS232/USB CONVERTER
T215	SQUARE PAN (not available for all models)
T216	SERIAL CABLE
T217	V-RANGE FUNCTION
T218	DENSITY FUNCTION'
T219	PAPER GRAMMAGE FUNCTION
T220	MAX LOAD DETERMINATION FUNTION
T221	TLP-50 PRINTER
T222	PROTECTION PLASTIC COVER

# 34 Warranty

- Duration of warranty is of 24 months from the date of purchase shown in the invoice concerning the product or by delivery note.
- Warranty covers all parts resulting defective at the origin. It does not cover mechanical or electronic parts damaged by wrong installation, tampering or incorrect use.
- Warranty does not cover damages caused by impacts, balance drops or drop of objects on weighing pan.
- Shipment to and from service centre is at customer charge.

# 35 Equipment disposal



This equipment is marked with the crossed out wheeled bin symbol to indicate that this equipment must not be disposed of with unsorted waste.

Instead it's your responsibility to correctly dispose of your equipment at lifecycle -end by handling it over to an authorized facility for separate collection and recycling. It's also your responsibility to decontaminate the equipment in case of biological, chemical and/or radiological contamination, so as to protect from health hazards the persons involved in the disposal and recycling of the equipment.

For more information about where you can drop off your waste of equipment, please contact your local dealer from whom you originally purchased this equipment.

By doing so, you will help to conserve natural and environmental resources and you will ensure that your equipment is recycled in a manner that protects human health.

Thank you

# Balances Series L / LG / LW / LGW

The models below indicated are all for internal use. Max altitude use: 4000m; Polltion grade: 2; Overvoltage category: II

Power sur	Power supply included.	. pc	INPLIT Switching	100-240Var~ F	INDLIT: Switching 100-240Vac 50/60Hz OLITPLIT:	1000 TOUM	ov DC 1000m4 Max absorbed nower 9//A	1 nower 9\\A					
Battery:			internal rechargea	ble (excluded i	internal rechargeable (excluded models with internal calibration)	I calibration)		5					
Enviromer adaption:	Enviroment condition adaption:	ſ	Selectable filters										
Autozero:			Selectable from Menu	enu									
Serial output:	out:		RS232C										
Operating range:	Operating temperature range:	re	+5°C - +35°C										
Model external version	Model internal calibration version	Weighing range(g)	(g) Resolution	(g) Tare ramge	Reproducib (standard (standard	(g) Linearity	Response time Baverage)	Span drift -10∴+30°C	Pan Dimension s (mm) s	Balance Dimension s (LxWxH) (mm)	Peso netto versione cal esterna (kg)	Net wight for internal calibration version (kg)	Minimum* Weight (class) for calibrating
L163 LG163	L163i LG163i	160	0.001	160	0.0006	<u>+</u> 0.003	≤ 3 sec.	±4 ppm/°C	110	345x215x155	3.5	3.5	g 100 (F1)
L203 LG203	L203i LG203i	210	0.001	210	0.0006	<u>+</u> 0.003	≤ 3 sec	±4 ppm/°C	110	345x215x155	3.5	3.5	g 100 (F1)
L303 LG303	L303i LG303i	310	0.001	310	0.0006	<u>+</u> 0.003	≤ 3 sec	$\pm 4 \text{ ppm/°C}$	110	345x215x155	3.5	3.5	g 200 (F1)
L423 LG423	L423i LG423i	420	0.001	420	0.001	<u>+</u> 0.004	≤4 sec	$\pm 4 \text{ ppm/°C}$	110	345x215x155	3.5	3.5	g 200 (F1)
L303D LG303D	L303Di LG303Di	150/310	0.001/0.01	150/310	0.0006/0.005	<u>+</u> 0.003/ <u>+</u> 0.02	≤ 3 sec	$\pm 4 \text{ ppm/°C}$	110	345x215x155	3.5	3.5	g 200 (F1)
LW163 LGW163	LW163i LGW163i	160	0.001	160	0.0006	<u>+</u> 0.003	≤ 3 sec.	±4 ppm/°C	110	345x215x345	5.5	5.5	g 100 (F1)

00

	Minimum* Weight (class) for calibrating	g 100 (F1)	g 200 (F1)	g 200 (F1)	g 200 (F1)	g 200 (F1)	g 200 (F1)	g 1000 (F1)	g 2000 (F1)	g 2000 (F1)	g 2000 (F1)	g 2000 (F1)	g 2000 (F1)
	Net wight for internal version (kg)	5.5	5.5	5.5	5.5	3.7	3.7	3.7	3.7	3.7			
	Peso netto versione cal.esterna (kg)	5.5	5.5	5.5	5.5	3.2	3.2	3.2	3.2	3.2	3.8	3.8	3.8
	Balance Dimension (HxWxH) s (mm)	345x215x345	345x215x345	345x215x345	345x215x345	345x215x100	345x215x100	345x215x100	345x215x100	345x215x100	345x215x110	345x215x110	345x215x110
	nsn9 Dimension Pan	110	110	110	110	160	160	160	160	160	160 (or 195x175)	160 (or 195x175)	195x175
	Span drift -10+30°C	±4 ppm/°C	±4 ppm/°C	±4 ppm/°C	±4 ppm/°C	±5 ppm/°C	±5 ppm/°C	±5 ppm/°C	±5 ppm/°C	±5 ppm/°C	±6 ppm/°C	±6 ppm/°C	±6 ppm/°C
	Response time average)	≤ 3 sec	≤ 3 sec	≤4 sec	≤ 3 sec	≤ 3 sec	≤ 3 sec	3 sec.	<u>&lt;</u> 3 sec.	$\leq$ 4 sec.	≤2 sec	≤2 sec	≤ 2 sec
	(g) Linearity	<u>+</u> 0.003	<u>+</u> 0.003	<u>+</u> 0.004	<u>+</u> 0.003/ <u>+</u> 0.02	± 0.02	± 0.02	± 0.03	± 0.03	± 0.04	± 0.1	± 0.1	± 0.1
c	Reproducil (standard (standard	0.0006	0.0006	0.001	0.0006/0.005	0.005	0.005	0.005	0.008	0.01	0.05	0.05	0.05
e	(g) Tare ramge	210	310	420	150/310	1000	1500	2200	3100	4200	4500	5200	6500
	(g) Kesolution	0.001	0.001	0.001	0.001/0.01	0.01	0.01	0.01	0.01	0.01	0.1	0.1	0.1
	Weighing range(g)	210	310	420	150/310	1000	1500	2200	3100	4200	4500	5200	6500
	Model internal calibration version	LW203i LGW203i	LW303i LGW303i	LW423i LGW423i	LW303Di LGW303Di	L1002i LG1002i	L1502i LG1502i	L2202i LG2202i	L3102i LG3102i	L4202i LG4202i	~	~	~
	Model external calibration version	LW203 LGW203	LW303 LGW303	LW423 LGW423	LW303D LGW303D	L1002 LG1002	L1502 LG1502	L2202 LG2202	L3102 LG3102	L4202 LG4202	L4501 LG4501	L5201 LG5201	L6501

Minimum* Weight (class) for calibrating		g 2000 (F1)	g 2000 (F1)	g 2000 (F1)	g 2000 (F1)	g 2000 (F1)
Peso netto versione cal.esterna (kg) internal calibration calibration version (kg)			3.8	3.8	3.8	3.8
Balance Dimension s (LxWxH) (mm)		345x215x110	345x215x110	345x215x110	345x215x110	345x215x110
Pan Dimension (mm) s		195x175	195x175	195x175	195x175	195x175
Span drift +10+30°C		± 6 ppm/°C	± 6 ppm/°C	± 6 ppm/°C	± 6 ppm/°C	± 6 ppm/°C
Response time Response		≤2 sec	≤ 2 sec.	≤ 3 sec.	≤ 3 sec.	≤4 sec.
(g) Linearity		± 0.2	<u>+</u> 0.2	<u>+</u> 0.2	<u>+</u> 0.3	<u>+</u> 0.4
Reproducib (standard (eviation)		0.05	0.05	0.08	0.08	0.1
(g) Tare ramge		8000	10000	12000	1 6000	20000
Resolution (g)		0.1	0.1	0.1	0.1	0.1
Weighing range(g)		8000	10000	12000	16000	20000
Model internal calibration version		/	/	/	/	~
Model external calibration version	LG6501	L8001 LG8001	L10001 LG10001	L12001 LG12001	L16001 LG16001	L20001 LG20001

\*Higher value weight can be used to calibrate the Balance by means of "LOAD" function (please refer to balance user manual)

List updated at: September 2011